



# Record video


When you selected all cameras, you see a preview in the main MediaRecorder window.

Image not found or type unknown

**TIP** If you selected multiple microphones, it can be useful to mute all but one audio channel in the preview. Select the audio source in the **Mute** column of the table under the preview. Muting audio in the preview does not influence the recording.

Image not found or type unknown


To start recording, press the **Start recording** button.

Image not found or type unknown

A pop-up window will appear, informing you about Noldus's Data Transparency Policy. Read it carefully before clicking the **I will inform the participants** button. You can also check the box **Don't show me again** to prevent further occurrence.

data transparency szek.png


The button then changes into the **Stop Recording** button and the Recording timer starts running.

Image not found or type unknown

Press the button again to stop recording. Recording stops automatically when you [specified a recording time](#) .

## how to access the recorded videos

To quickly access the recorded videos, press the **Open target folder** button in the main MediaRecorder window.

Image not found or type unknown

The video files are stored in the default video file location, or the location you specified. See [To select the video file location](#).