



Record video

When you selected all cameras, you see a preview in the main MediaRecorder window.

Image not found or type unknown

TIP If you selected multiple microphones, it can be useful to mute all but one audio channel in the preview. Select the audio source in the **Mute** column of the table under the preview. Muting audio in the preview does not influence the recording.

Image not found or type unknown


To start recording, press the **Start recording** button.

Image not found or type unknown

A pop-up window will appear, informing you about Noldus's Data Transparency Policy. Read it carefully before clicking the **I will inform the participants** button. You can also check the box **Don't show me again** to prevent further occurrence.

data transparency szek.png

The button then changes into the **Stop Recording** button and the Recording timer starts running.

Image not found or type unknown

Press the button again to stop recording. Recording stops automatically when you [specified a recording time](#) .

how to access the recorded videos

To quickly access the recorded videos, press the **Open target folder** button in the main MediaRecorder window.

Image not found or type unknown

The video files are stored in the default video file location, or the location you specified. See [To select the video file location](#).