

Change video order in preview

Aim

To change the order of the video cameras, for example to change the position of the images in a Picture in Picture recording.

Prerequisite

You selected the recording devices in the **Video Settings** window.

See [Select a video device](#)

Procedure

1. Drag and drop rows. in the table with video devices or drag and drop the preview images.
2. Click the **Refresh** button.

See [Refresh the preview](#)

Note

You cannot drag and drop Picture in Picture, or Picture by Picture preview images. Drag and drop the rows in the table with video devices instead.

Revision #5

Created 29 November 2023 21:40:09

Updated 20 December 2023 12:47:08