

# Change video order in preview

## Aim

To change the order of the video cameras, for example to change the position of the images in a Picture in Picture recording.

## Prerequisite

You selected the recording devices in the **Video Settings** window.

See [Select a video device](#)

## Procedure

1. Drag and drop rows. in the table with video devices or drag and drop the preview images.
2. Click the **Refresh** button.

See [Refresh the preview](#)

## Note

You cannot drag and drop Picture in Picture, or Picture by Picture preview images. Drag and drop the rows in the table with video devices instead.

---

Revision #5

Created 29 November 2023 22:40:09

Updated 20 December 2023 13:47:08