



# Change video standard

## Aim

To specify the correct video standard to be used by the Picolo Alert PCIe card.

## Procedure

1. Open **Video settings** the first time you select the Picolo Alert PCIe card.  
nd or type unknown
2. Select the correct video standard, depending on your cameras. It's typically PAL for Europe and NTSC for the US. You need to do this four times, selecting all four channels from the **Board** drop-down box (channels are numbered #0-#3). Press **Apply** after setting the correct Video Standard for each channel. nd or type unknown

You should see the preview after setting the Video Standard correctly. Failing this will result in a black preview. You can only modify this setting the first time before the preview is running, after that it will become disabled. To be able to modify it, simply deselect the camera from the video inputs, and then select it again, to make the setting editable once more.

---

Revision #3

Created 29 November 2023 23:52:43

Updated 2 January 2024 17:51:01