

Configure the Basler USB 3 camera

Aim

To adjust settings in the USB 3 camera, like the video frame rate, resolution, and exposure.

NOTE The camera settings are stored by the camera itself, not on the PC. When you work with multiple cameras, repeat the procedure for each camera.


Procedure

1. Make sure MediaRecorder (and EthoVision XT, when present on the computer) are not running.
2. Start the **Basler Pylon Viewer** software that comes with the USB 3 camera.

 Image not found or type unknown

NOTE The Basler Pylon Viewer software is installed automatically when you install the camera drivers.

3. In the **Devices** panel, under **USB**, double-click the Basler camera.

 Image not found or type unknown

4. Under **Features**, open the item **Basler acA1920-155um** and choose the settings you would like to change. For details, see the corresponding settings in [Configure the cameras](#) in [Set up GigE cameras](#).
5. Save the settings under **User Set Control**.

See also

- [Supported USB cameras](#) in [Recording devices](#)

Revision #7

Created 1 December 2023 11:21:38

Updated 2 January 2024 21:45:42