

# Configure the Basler USB 3 camera

## Aim

To adjust settings in the USB 3 camera, like the video frame rate, resolution, and exposure.

**NOTE** The camera settings are stored by the camera itself, not on the PC. When you work with multiple cameras, repeat the procedure for each camera.


## Procedure

1. Make sure MediaRecorder (and EthoVision XT, when present on the computer) are not running.
2. Start the **Basler Pylon Viewer** software that comes with the USB 3 camera.

 Image not found or type unknown

**NOTE** The Basler Pylon Viewer software is installed automatically when you install the camera drivers.

3. In the **Devices** panel, under **USB**, double-click the Basler camera.

 Image not found or type unknown

4. Under **Features**, open the item **Basler acA1920-155um** and choose the settings you would like to change. For details, see the corresponding settings in [Configure the cameras](#) in [Set up GigE cameras](#).

5. Save the settings under **User Set Control**.

## See also

- [Supported USB cameras](#) in [Recording devices](#)

---

Revision #7

Created 1 December 2023 10:21:38

Updated 2 January 2024 20:45:42