

# Ethernet card setup

## Aim

To insert the 1 Gb Ethernet card into your computer and to configure it to work with MediaRecorder.

[image.png](#) image not found or type unknown

## Prerequisites

- You have IP or GigE cameras.
- You have an Intel PRO/1000 GT or Intel Pro/ 1000 PT Ethernet card.
- You purchased the computer with MediaRecorder yourself, or you bought the cameras after you received the computer.

If you bought a complete system from Noldus IT, the card has been installed and set up. You can skip this topic.

## Procedure

Carry out the following steps:

1. [Install the Ethernet card](#)
2. [Set Ethernet card IP address](#)

---

Revision #6

Created 1 December 2023 09:31:28

Updated 2 January 2024 18:14:34