

# Install the drivers for the Picolo Alert card

## Aim

To install the software that allows communication between the Euresys card and the computer operating systems.

## Prerequisites

- You use MediaRecorder with analog cameras.
- You did not buy a complete setup from Noldus IT.
- You installed the Euresys Picolo Alert card into the computer.

## Procedure

1. Insert the MediaRecorder installation USB stick into your computer.
2. If you have the Picolo Alert card, open the folder Drivers and Tools\Drivers\Euresys\Alert and double-click the file **multicam-win10-6.15.1.3573.exe**
3. Follow the instructions on your screen.
4. When installation is finished, a message appears that the drivers were successfully installed. Click **Close**.
5. Installing the Picolo card drivers involves a a computer restart. After restart, the installation is finished.

Continue with installing MediaRecorder. See [Installation](#)

---

Revision #5

Created 29 November 2023 22:50:53

Updated 20 December 2023 12:47:08