

# Record your videos

## Aim

To create video files from the video stream of the connected recording devices.

## Prerequisites

- Your devices are connected to the computer with MediaRecorder, and, if applicable, switched on.
- MediaRecorder is open.
- You selected the cameras and, optionally, microphones and specified the output of the video files in the **Video Settings** window (**File > Video Settings**).

## Procedure

1. When you selected all settings in the **Video Settings** window, click **OK** to return to the Main window. You will see a preview of the cameras.

 image.png and or type unknown

2. Optionally, adjust options like the video file location, the video file base name or the order of the video files. See [Options](#)
3. Press the **Start recording** button.

 image.png and or type unknown

4. A pop-up window will appear, informing you about Noldus's Data Transparency Policy. Read it carefully before clicking the **I will inform the participants** button. You can also check the box **Don't show me again** to prevent further occurrence.

 data transparency szerk.png and or type unknown

5. The button changes into the **Stop Recording** button and the Recording timer starts running.

 image.png and or type unknown

6. Press the button again to stop recording. Recording stops automatically when you specified a [Maximum recording time](#).

## Note

If you have IP cameras, see [Pan, Tilt, Zoom control with IP cameras](#) how to change the camera view before or during recording.