

Set video file base name

Aim

To select a word or phrase to start the video file name with.

The video file name consists of the **File base** name followed by:

- The video input name.
- The date and the time the videos were recorded.
- The number of the video camera.

Example – The **File base name** is Video, you changed the name of Video 2 to Webcam in the **Video Settings** window (**File > Video Settings**), the date is 11 May 2016, the time 10:54:43 and you create 2 videos simultaneously. The name of the second video will then be:

Video Webcam 5_11_2016 10_54_43 AM 2.mp4

Procedure

Enter a name in the **File base name** field.

Image not found or type unknown

Note

The date and time depend on your computer's regional settings. To change these, open the **Control panel** and then **Region and Language**. Select your preferences in one of the tabs.

Revision #3

Created 29 November 2023 21:37:59

Updated 20 December 2023 23:01:20