

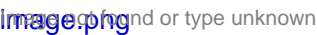
Show frame rate, resolution and format

Aim

To display hidden rows in the table with video devices.

Procedure

Right-click one of the column headers and select the preferred options.

Image not found or type unknown

Notes

- The **Format** column shows the color space of the video file. The column is hidden by default, because the optimal color space is selected by default. See the following page for more information on color space.
<http://www.noldus.com/knowledge-base/what-color-space-format-used-media-recorder>.
- **TIP** Follow the same procedure to show or hide the Frame rate, Resolution, and Format columns in the **Video Settings** window.