

Synchronize videos with events

Aim

To synchronize the MediaRecorder videos with the event log of The Observer XT.

When is manual synchronization not needed?

Videos created by MediaRecorder are automatically synchronized with the events in The Observer XT when you followed the procedure in [MediaRecorder with The Observer XT](#).

An automatic offset is applied to the videos. See [View the automatic offset of videos imported into The Observer XT](#) to see which offset was applied.

When is manual synchronization needed?

You need to synchronize videos and events manually in the following situations:

- You start video recording manually and import the videos into The Observer XT.
- You have an older version of The Observer XT that does not support N-Linx data transfer.
- See [Synchronize manually](#) for the procedure.

Revision #5

Created 29 November 2023 23:25:08

Updated 20 December 2023 13:47:09