

# The steps to install MediaRecorder

## IMPORTANT

**If you have an older version of MediaRecorder** - Uninstall MediaRecorder 6.0 and older versions before you install MediaRecorder 6.5.

**If you have a hardware key** - Do not insert your key before the installation is complete. If you do so, other drivers are automatically installed, which may lead to incorrect functioning of the license key. Drivers of the license key should be installed using the installation USB stick.

See [Install MediaRecorder and device drivers](#)

## Procedure

1. [Turn off automatic updates for device drivers](#)
2. [Select correct power options](#)
3. [Install MediaRecorder and device drivers](#)

---

Revision #8

Created 29 November 2023 17:18:13

Updated 20 December 2023 22:21:58