

# The steps to install USB 3 cameras

## Prerequisite

You have USB 3 cameras that supported with MediaRecorder.

See [Supported USB cameras](#) in [Recording devices](#)

## Which situation applies to you?

- You ordered the computer, cameras and MediaRecorder from Noldus IT.

In case of Basler USB 3.0 cameras, your computer has a USB 3 card and its drivers are installed. The cameras are configured. Also MediaRecorder has been set up to use with these cameras. You can connect your cameras and start recording.

See [Connect the USB 3 cameras to the PC](#) and [Record video](#)

- You bought the computer somewhere else.

Carry out the entire [Procedure](#) below.

## Procedure

1. In case of Basler USB 3 cameras, install the supported USB 3 interface card, even if your computer already has USB 3 ports.

See [Install the USB 3 interface card](#)

2. Carry out the prerequisite steps before installing MediaRecorder.

See [The steps to install MediaRecorder](#) in [Installation](#)

3. Install (MediaRecorder with) the camera driver. Installing the camera driver is only needed for the Basler USB 3 camera. For the Logitech BRIO no driver installation is needed.

See [Install the USB 3 camera driver](#)

4. [Connect the USB 3 cameras to the PC](#)
5. [Configure the Basler USB 3 camera](#)

## See also

[Supported USB cameras](#) in [Recording devices](#)