

# Upgrade MediaRecorder

## Aim

To upgrade an older version of the MediaRecorder software to version 6.5.

## Prerequisites

You have an older MediaRecorder version.

You obtained an activation key for your MediaRecorder license.

## Procedure

1. Download the installation files from [my.noldus.com](https://my.noldus.com), and extract the contents.
2. Browse to the file **MediaRecorder [Version number] Setup.exe** and double-click it.
3. Follow the steps on your screen to install MediaRecorder 6.5.
4. [Activate your MediaRecorder license](#).

---

Revision #9

Created 29 November 2023 17:07:09

Updated 21 December 2023 10:20:34 by Niek Wilmink