

Customize the Session window

What do you want to do?

- [Dock/Undock Event log or Time Line](#)
- [Dock/Undock panes and tabs](#)
- [Display a selection of events](#)
- [Collapse all subjects](#)

Dock/Undock Event log or Time Line

You can undock and move the Event log and Time Line window, for example to display them on a separate monitor. To dock the windows again, drag the window to the main Viso window. Docking buttons appear. Move the window to one of these docking items. A blue rectangle now shows where the window will be positioned. Release the mouse button.

[image.png](#) and or type unknown

Dock/Undock panes and tabs

You can undock the pane with the **Events**, **Codes** and **Participant** tabs, and the individual tabs. It can, for example, be useful to move the **Codes** tab to a separate monitor.

Docking/undocking of panes and tabs applies to Viso Viewer and is implemented for all users.

To undock a pane with its tabs

Click the title and drag it to a new position.

[image.png](#) and or type unknown

To dock a pane again, drag it to the main Viso window. Docking buttons appear. Move the pane to one of these docking items. A blue rectangle now shows where the pane will be positioned. Release the mouse button.

[image.png](#) and or type unknown

To undock a separate tab

Click the tab at the bottom of the pane and drag it to a new position.

[image.png](#) and or type unknown

To dock the tab again, drag it to the main Viso window. Two types of docking buttons appear.

If you release the mouse button while it is positioned on the docking button shown in the figure below, it is redocked in the pane.

[image.png](#) and or type unknown

If you release the mouse button while it is positioned on the docking button shown in the figure below, the separate tab is docked in the Viso window.




[image.png](#) and or type unknown

Display a selection of events

aim


To select a subset of events in the Event log and, optionally, in the Timeline. This can be useful to, for example, quickly display all behaviors of one subject, or a specific behavior of all subjects.

procedure

1. Open the **Events** pane.
2. Click the filter button next to **Subject** or **Marker**.

3. Specify the selection. The result is shown in the Event log.

The filter is also applied in the Timeline.


To disable the filter in the Timeline

To see in what context the selected events took place, you may want to display all events in the Timeline, instead of only the ones specified in the Event log filter. To do so:

1. Right-click the Timeline.
2. Select **Show all events**.


NOTE To restore the filter in the Timeline, right-click it and select **Show filtered events only**.

Collapse all subjects

aim

To quickly collapse all subjects in the Timeline, so that no markers are shown. You can then expand the subjects that you are interested in to view only their markers.



procedure

Right-click the Timeline and select **Collapse all subjects**.

NOTE To expand all collapsed subject again, right-click the Timeline and select **Expand all subjects**.

Revision #4

Created 15 January 2024 11:38:50

Updated 11 March 2024 14:26:38