

# Export annotations to text file

## aim

To export the annotations to a Comma Separated Value text file (CSV-file) so that you can analyze them in another application, like Excel.

## prerequisites

- Your license includes the Annotation Module.
- You created the session, have received session ownership or are a Technician or Administrator.

## procedure

1. Open the **Sessions** tab.
2. Select the sessions to export.
3. Right-click one of the sessions and select **Export to CSV**.
4. Choose a location and click **Export**.

**NOTE** The export file is named after the imported session, with the location name, and the date and time the session was created added to the filename.

---

Revision #2

Created 15 January 2024 10:38:50

Updated 10 March 2024 18:38:40