


# How Viso annotations appear in The Observer XT


## In the Coding Scheme

In The Observer XT, choose **Setup > Open Coding Scheme**, to open the Coding Scheme. Viso subjects appear in the **Subjects** panel. Viso markers are added as Point events in a Start-Stop group. Remarks are not added to the Coding Scheme.

 Image not found or type unknown

## In the observation

Choose **Observe > Observation > Open > [session name and time]** to view the annotations in the Event Log. Viso Subjects are also Subjects in The Observer XT. Markers are point events and are indicated with blue dots. Remarks are imported as comments in the Event Log.

 Image not found or type unknown

To jump to the video frame at which a certain marker or remark was scored, click its row in the Event Log.

---

Revision #1

Created 15 January 2024 10:38:52

Updated 15 January 2024 10:43:02