

# Reclaim the role of annotator


## aim

To claim back the role of annotator that you gave to another user, for instance, a colleague or one of your students.

## prerequisites

- Your license includes the Annotation Module.
- You created the session or you received session ownership or you are a Technician or Administrator.

## procedure

1. Open the **Sessions** tab to view the list of all the sessions.
2. Right-click the appropriate session and select **Change annotator**.  
The screenshot shows a dialog box with the title 'Change annotator'. It contains a text input field with the placeholder text 'Image not found or type unknown'. Below the input field is a button labeled 'OK'.
3. Select your name and click **OK**.

---

Revision #2

Created 15 January 2024 11:38:50

Updated 8 March 2024 16:37:56