

Reclaim the role of annotator

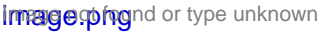
aim

To claim back the role of annotator that you gave to another user, for instance, a colleague or one of your students.

prerequisites

- Your license includes the Annotation Module.
- You created the session or you received session ownership or you are a Technician or Administrator.

procedure

1. Open the **Sessions** tab to view the list of all the sessions.
2. Right-click the appropriate session and select **Change annotator**.
 Image copied and type unknown
3. Select your name and click **OK**.

Revision #2

Created 15 January 2024 11:38:50

Updated 8 March 2024 16:37:56